

# UltraMarathonRT<sup>®</sup>

20,000 U/mL | Store at -20°C

## TWO-STEP RT-qPCR KIT

This kit uses a green dye that is compatible with most of the real-time PCR systems available (Ex 498/Em 522), providing sensitive detection for any difficult RNA templates regardless of PCR amplicon location.

### Components Provided

UltraMarathonRT

2x RT Reaction Buffer Q

dNTP Mix

Oligo (dT)<sub>18</sub> Primer

Random Primer (15 mer)

2x qPCR Green Dye Master Mix

Nuclease-Free Water



For product details, visit: [www.RNAConnect.com](http://www.RNAConnect.com)

# RT-qPCR Quick Start Protocol

---

## ANNEALING RT PRIMERS TO RNA TEMPLATES

1. Gently mix the following components in a nuclease-free microcentrifuge tube by tapping the tube and collect the liquid with a quick spin.

Components	Volume (6 $\mu$ L total)
0.1 pg - 2 $\mu$ g of total cellular RNA or 0.1 pg - 500 ng of mRNA	'x' $\mu$ L
Oligo (dT) <sub>18</sub> (5 $\mu$ M), or Randomer (15 mer) (10 $\mu$ M), or Gene-specific primer (2 $\mu$ M)	1 $\mu$ L
dNTP Mix (10 mM each)	1 $\mu$ L
Nuclease-Free Water	(4 - 'x') $\mu$ L

---

2. Incubate at 95°C for 30 sec and then snap cool on ice to anneal the primer to the template.

## PREPARING THE RT REACTION MIXTURE

3. Gently mix the following components in a nuclease-free microcentrifuge tube by tapping the tube and collect the liquid with a quick spin.

Components	Volume (14 $\mu$ L total)
2x RT Reaction Buffer Q	10 $\mu$ L
UltraMarathonRT	1 $\mu$ L
RNase Out™ (40 U/ $\mu$ L) (optional)	1 $\mu$ L
Nuclease-Free Water	2 $\mu$ L

---

4. Add RT reaction mix to the annealed RNA (step 2) and mix gently by tapping the tube.
5. Incubate at 30°C for 15 min to carry out reverse transcription. For RNA > 12 kb, 30°C for 20 - 60 min is recommended.
6. Inactivate the enzyme by heating at 95°C for 1 min.
7. The cDNA can be stored at -20°C or immediately used for PCR amplification.

## AMPLIFY THE cDNA USING qPCR

8. Gently mix the following components to create a single qPCR reaction.

Components	Volume (20 $\mu$ L total)
2x qPCR Green Dye Master Mix	10 $\mu$ L
Forward primer (10 $\mu$ M)	0.6 $\mu$ L
Reverse primer (10 $\mu$ M)	0.6 $\mu$ L
Template DNA	1 - 2 $\mu$ L of uMRT cDNA (step 7)
PCR grade water	Added to 20 $\mu$ L

→ Up to 2  $\mu$ L of the unpurified cDNA product (step 7) can be used for a 20  $\mu$ L qPCR reaction. More than 2  $\mu$ L of the unpurified cDNA may inhibit the PCR reaction.

### qPCR DATA COLLECTION:

Detection	Excitation Peak	Emission Peak
Green dye reagents	498 nm	522 nm

→ Performing a melting curve analysis is important to evaluate the specificity of any dye-based qPCR assay. Specificity can also be confirmed using electrophoresis analysis on an agarose gel.

## qPCR CYCLING CONDITIONS:

Stage of Cycle		Temp (°C)	Time
Pre-Denaturation		98	1:30 min
Amplification (40-45 cycles)	Melt	98	10 sec
	Anneal	60	10 sec
	Extension <sup>1</sup>	68	30 sec
Melting Curve	Step 1	95	5 sec
	Step 2	65	1:00 min
	Step 3 <sup>2</sup>	97	-
Cooling		40	10 sec

1. Be sure acquisition is set to 'single' at the Extension step. An extension time of 30 seconds is sufficient for all amplicon sizes  $\leq 500$  bp.
  2. Be sure acquisition is set to 'continuous' with a ramp rate of  $0.05$  °C / sec at Step 3 of the Melting Curve cycle for smooth curves. Ramp rate for all other stages can be set to their maximum value.
- For detailed qPCR troubleshooting suggestions, please view our full protocol available online (see QR code on cover).

## STORAGE CONDITIONS

uMRT is stable in  $-20^{\circ}\text{C}$  for up to 3 months. For long-term storage,  $-80^{\circ}\text{C}$  is recommended.

This product and its labels and packaging may be covered by one or more patents, trademarks or copyrights owned or licensed by RNAConnect, Inc. The use of this product may require additional third-party intellectual property rights for certain applications. This product is for research use only and is not intended for use in humans. All product names and trademarks are the property of their respective owners. © 2025 RNAConnect, Inc. All rights reserved.